

HLA Positive Control

(Rabbit Anti-Human Leucocyte Serum)

Package Insert

PC-HLA 0.5 ml
SPC-HLA 250 µl

INTENDED USE

For use as a control that reacts with human leukocytes.

For Research Use only.

SUMMARY/EXPLANATION

This HLA Positive Control is an anti-human leukocyte serum prepared from rabbits that have been immunized with human thymocytes.

- Store frozen at $\leq -65^{\circ}\text{C}$.
- Repeated thawing and freezing is detrimental to antibody.

Warning: Animal blood products have the potential to transmit diseases. Take appropriate precautions.

Caution: This product contains <0.1% sodium azide. Sodium azide reacts with lead and copper plumbing forming highly explosive metal azides.

Dispose of according to local regulations.

INSTRUMENTS/EQUIPMENT

(Not included with reagent)

$\leq -65^{\circ}\text{C}$ Freezer

INSTRUCTIONS FOR USE

1. Thaw prior to use.
2. Aliquot into microcentrifuge tubes in volume suitable for single use.
3. Freeze aliquots at $\leq -65^{\circ}\text{C}$.
4. Thaw only as needed for use.

REACTIVITY

Determine reactivity in your laboratory's test system.

STORAGE, HANDLING, SHIPPING

HLA Positive Control is shipped frozen, on dry ice. Promptly store the frozen serum at $\leq -65^{\circ}\text{C}$, upon arrival.

HLA Positive Control
Rabbit Anti-Human Leucocyte Serum



GTi DIAGNOSTICS®
Good science starts with people.®

- FOR RESEARCH USE ONLY
- STORE AT $\leq -65^{\circ}\text{C}$

20925 Crossroads Circle, Suite 200
Waukesha, WI 53186-4054
(262) 754-1000 or 1-800-233-1843

Cat. NO.: PC-HLA, SPC-HLA
Rev. 12 October 2007

www.gtidiagnostics.com